

SYSTEC CONNECT SAVE TO FOLDER

for Systemec DX-Series, Systemec VX-Series and Systemec HX-Series (2D) Autoclaves

SAVE TO FOLDER OPTION

Automatic data transfer via FTP/SFTP/FTPS. Process data is automatically saved to a user-definable folder in the computer network (e.g. to a server, to a PC or to other storage devices connected to the network). Process data error logs and Audit Trail can be exported manually to the same folder in the computer network.



PARAMETERS

FTP-Export

Parameter	FtpCSV
Description	Enables the possibility to export CSV files of the process log, error log and Audit Trails to an FTP server.
Unit	Minimum Maximum Default
-	- - off

Parameter	FtpPDF
Description	Enables the possibility to export PDF files of the process log (without tabular listing of the measured values), error log and audit trails to an FTP server.
Unit	Minimum Maximum Default
-	- - off

Parameter	FtpExtPDF
Description	Enables the possibility to export extended PDF files of the process log including sensor statistics, program parameters and tabular listing of the measured values to a FTP server.
Unit	Minimum Maximum Default
-	- - off

Parameter	FtpURL
Description	Defines the address of the FTP server to which the PDF/CSV documentation should be exported.
Unit	Minimum Maximum Default
Text	- - -

Parameter	FtpUser
Description	Defines the username for the FTP server.
Unit	Minimum Maximum Default
Text	- - -

Parameter	FtpPW
Description	Defines the password for the FTP server.
Unit	Minimum Maximum Default
Text	- - -

PRICE & AVAILABILITY

Product name	Product description	Art. No.	Price*	Availability
Save to Folder	Automatic data transfer via FTP/SFTP/FTPS	036-1	see price list	now

Subject to change and errors excepted.
*excl. VAT.